

# iDTECT® INTEGRATION SITE ANALYSIS CLONE SELECTION AND GENETIC STABILITY WITH NEXT GENERATION SEQUENCING (NGS): COMPLIANT, SAFER AND FASTER.

## MODALITIES

- **Engineered Cell Lines/Cell Therapies:**
  - CHO/mAbs
  - Human cells/lentivirus
  - Engineered T-cells
- **Cultured meat cell banks**

## WHY CHOOSE iDTECT® INTEGRATION SITE ANALYSIS

- ✓ Better understanding of your clones
- ✓ Resolve concatemers, truncations and inverted repeats more easily than short-read integration site analysis

## COMPARISON OF INTEGRATION SITE ANALYSIS METHODS

	 <b>FISH</b>	 <b>Sanger Sequencing</b>	 <b>Short-read NGS</b>	 <b>Long-read ISA</b>
<b>Integration Site Resolution</b>	☆☆☆☆☆	★★★★☆	★★★★★	★★★★★
<b>Transgene Location</b>	★★★★☆	★★★★☆	★★★★★	★★★★★
<b>Concatemer &amp; Inverted Repeats</b>	☆☆☆☆☆	☆☆☆☆☆	★★★★☆	★★★★★

As part of the development of a cell line for manufacturing, candidate clones are selected based on a number of attributes including production yield, number of transgene copies, as well as the structure and location of the transgene in the host cell genome. Indeed, a certain level of genetic characterization of the cell line is expected by the regulatory authorities. As such, developers can benefit from a deeper understanding of the integration site of the transgene when making clone selection. For example, transgenes located in areas of instability or regions that are prone to epigenetic silencing may lead to a significant reduction in manufacturing yields over the cell line's manufacturing lifetime.

Classically, integration site analysis has relied on low resolution techniques such as fluorescent *in situ* hybridization (FISH) or southern blotting. Higher resolution approaches such as Sanger sequencing provide the required precision, but can struggle to identify complex integration events. A targeted analysis based on short-

read NGS such as Targeted Locus Amplification (TLA) uses genome crosslinking to generate fragments up to 100kb either side of a known primer site. This method provides some information on insertion site, but still struggles with identification of concatemers, truncations and inverted repeats.

To provide increased resolution, PathoQuest has developed a targeted, nanopore sequencing driven approach for Integration Site Analysis (ISA). This NGS assay based on Oxford Nanopore's long-read sequencing technology, further enhances the accuracy and reliability of genetic characterization. By targeting regions within the transgene expression cassette, both the transgene and it(s) associated junction site(s) can be comprehensively sequenced. Long read nanopore sequencing has the advantage over short-read ISA methods as it better characterizes concatemers, truncations and inverted repeats, providing a superior and more accurate result.

SAMPLE REQUIREMENTS	SHIPMENT & STORAGE	TURNAROUND TIME (CALENDAR DAYS)	DELIVERABLES TO CLIENT
5x10 <sup>6</sup> cells (≥10µg DNA per sample) Backup sample required BSL1 or 2*	Dry Ice / -80°C	· Standard: 35 days** · Fast Track: contact us	· Report · Raw Data

\* Biosafety level classifications can vary between regulatory authorities – contact PathoQuest to discuss.  
\*\* Turnaround time per round.



To find out how Pathoquest, a Charles River Company can better characterize your cells, visit:  
[www.pathoquest.com/services/integration-site-analysis](http://www.pathoquest.com/services/integration-site-analysis)

BioPark – BatB 11 Rue  
Watt 75013, Paris

466 Devon Park Dr,  
Wayne, PA 19087, USA